

Density Standards - Premium Range

Summary of Features & Benefits:

Commercial Benefits

- Extensive range (0.6407 - 3.1140)
- No toxic heavy metals used in any formulation
- Can be used with any brand or type of density measuring instrument
- Presented in a high quality tamper proof amber glass bottle
- Customised formulations available

Technical Benefits

- High accuracy products
- Tested using a fundamental measurement technique (Bingham Pycnometer)
- All products tested in accordance with ASTM D1480 Guidelines
- Consistency of product - Independent, Traceable, Certified
- Certificates of Analysis and Safety Data Sheets available online

Reagecon manufactures an extensive range of Density Standards in accordance with ASTM D1480-12 for testing of Density or Relative Density (specific and API gravity) by Bingham Pycnometer. These materials can be used as calibration standards for density measurement by pycnometric techniques, vibrational techniques or hydrometer based techniques.

The product range includes:

- 0.6960 - 3.1140g/ml @ 15°C
- 0.6619 - 3.1096g/ml @ 20°C
- 0.6878 - 3.1043g/ml @ 25°C
- 0.6752 - 3.0852g/ml @ 40°C
- 0.6668 - 3.0721g/ml @ 50°C
- 0.6582 - 1.0478g/ml @ 60°C
- 0.6407 - 1.0302g/ml @ 80°C

The products are prepared gravimetrically on a weight/weight basis. Both solute and solvent are weighed on a balance calibrated by Reagecon engineers using OIML traceable weights. Reagecon holds ISO 17025 accreditation for calibration of laboratory balances (INAB Ref:265C). The resulting Balance Certificate of Calibration is issued in accordance with the requirements of ISO/IEC 17025. The density of each standard is established using a high performance set of fully calibrated reference pycnometers.



Density Standards @ 15°C

Product No.	Description	Pack Size
DEN15010PY	Density Standard 0.6960g/ml @15°C	100ml
DEN15020PY	Density Standard 0.7073g/ml @15°C	100ml
DEN15030 PY	Density Standard 0.7184g/ml @15°C	100ml
DEN15040PY	Density Standard 0.7298g/ml @15°C	100ml
DEN15050PY	Density Standard 0.7411g/ml @15°C	100ml
DEN15060PY	Density Standard 0.7524g/ml @15°C	100ml
DEN15070PY	Density Standard 0.7721g/ml @15°C	100ml
DEN15080PY	Density Standard 0.7933g/ml @15°C	100ml
DEN15090PY	Density Standard 0.8168g/ml @15°C	100ml
DEN15100PY	Density Standard 0.8428g/ml @15°C	100ml
DEN15110PY	Density Standard 0.8715g/ml @15°C	100ml
DEN15120PY	Density Standard 0.9135g/ml @15°C	100ml
DEN15130PY	Density Standard 0.9514g/ml @15°C	100ml
DEN15140PY	Density Standard 1.0040g/ml @15°C	100ml
DEN15150PY	Density Standard 1.0337g/ml @15°C	100ml
DEN15160PY	Density Standard 1.0828g/ml @15°C	100ml
DEN15170PY	Density Standard 1.1661g/ml @15°C	100ml
DEN15180PY	Density Standard 1.2498g/ml @15°C	100ml
DEN15190PY	Density Standard 1.3318g/ml @15°C	100ml
DEN15200PY	Density Standard 1.4152g/ml @15°C	100ml
DEN15210PY	Density Standard 1.5820g/ml @15°C	100ml
DEN15215PY	Density Standard 1.6459g/ml @15°C	100ml
DEN15220PY	Density Standard 1.7495g/ml @15°C	100ml
DEN15225PY	Density Standard 1.8366g/ml @15°C	100ml
DEN15230PY	Density Standard 1.9171g/ml @15°C	100ml
DEN15240PY	Density Standard 2.0846g/ml @15°C	100ml
DEN15250PY	Density Standard 2.2568g/ml @15°C	100ml
DEN15260PY	Density Standard 2.4261g/ml @15°C	100ml
DEN15270PY	Density Standard 2.6055g/ml @15°C	100ml
DEN15280PY	Density Standard 2.7588g/ml @15°C	100ml
DEN15290PY	Density Standard 2.9418g/ml @15°C	100ml
DEN15300PY	Density Standard 3.1140g/ml @15°C	100ml

Density Standards @ 20°C

Product No.	Description	Pack Size
DEN20010PY	Density Standard 0.6919g/ml @20°C	100ml
DEN20020PY	Density Standard 0.7033g/ml @20°C	100ml
DEN20030PY	Density Standard 0.7148g/ml @20°C	100ml
DEN20040PY	Density Standard 0.7261g/ml @20°C	100ml
DEN20050PY	Density Standard 0.7374g/ml @20°C	100ml
DEN20060PY	Density Standard 0.7488g/ml @20°C	100ml
DEN20070PY	Density Standard 0.7683g/ml @20°C	100ml
DEN20080PY	Density Standard 0.7893g/ml @20°C	100ml
DEN20090PY	Density Standard 0.8126g/ml @20°C	100ml
DEN20100PY	Density Standard 0.8384g/ml @20°C	100ml
DEN20110PY	Density Standard 0.8668g/ml @20°C	100ml
DEN20120PY	Density Standard 0.9098g/ml @20°C	100ml
DEN20130PY	Density Standard 0.9476g/ml @20°C	100ml
DEN20140PY	Density Standard 1.0005g/ml @20°C	100ml
DEN20150PY	Density Standard 1.0301g/ml @20°C	100ml
DEN20160PY	Density Standard 1.0792g/ml @20°C	100ml
DEN20170PY	Density Standard 1.1651g/ml @20°C	100ml
DEN20180PY	Density Standard 1.2486g/ml @20°C	100ml
DEN20190PY	Density Standard 1.3304g/ml @20°C	100ml
DEN20200PY	Density Standard 1.4136g/ml @20°C	100ml
DEN20210PY	Density Standard 1.5799g/ml @20°C	100ml
DEN20220PY	Density Standard 1.7470g/ml @20°C	100ml
DEN20230PY	Density Standard 1.9141g/ml @20°C	100ml
DEN20240PY	Density Standard 2.0812g/ml @20°C	100ml
DEN20250PY	Density Standard 2.2531g/ml @20°C	100ml
DEN20260PY	Density Standard 2.4219g/ml @20°C	100ml
DEN20270PY	Density Standard 2.6011g/ml @20°C	100ml
DEN20280PY	Density Standard 2.7542g/ml @20°C	100ml
DEN20290PY	Density Standard 2.9370g/ml @20°C	100ml
DEN20300PY	Density Standard 3.1096g/ml @20°C	100ml

Density Standards @ 25°C

Product No.	Description	Pack Size
DEN25010PY	Density Standard 0.6878g/ml @25°C	100ml
DEN25020PY	Density Standard 0.6993g/ml @25°C	100ml
DEN25030PY	Density Standard 0.7111g/ml @25°C	100ml
DEN25040PY	Density Standard 0.7223g/ml @25°C	100ml
DEN25050PY	Density Standard 0.7337g/ml @25°C	100ml
DEN25060PY	Density Standard 0.7452g/ml @25°C	100ml
DEN25070PY	Density Standard 0.7645g/ml @25°C	100ml
DEN25080PY	Density Standard 0.7853g/ml @25°C	100ml
DEN25090PY	Density Standard 0.8084g/ml @25°C	100ml
DEN25100PY	Density Standard 0.8340g/ml @25°C	100ml
DEN25110PY	Density Standard 0.8622g/ml @25°C	100ml
DEN25120PY	Density Standard 0.9060g/ml @25°C	100ml
DEN25130PY	Density Standard 0.9438g/ml @25°C	100ml
DEN25140PY	Density Standard 0.9969g/ml @25°C	100ml
DEN25150PY	Density Standard 1.0265g/ml @25°C	100ml
DEN25160PY	Density Standard 1.0755g/ml @25°C	100ml
DEN25170PY	Density Standard 1.1639g/ml @25°C	100ml
DEN2512PY	Density Standard 1.2000g/ml @25°C	100ml
DEN25180PY	Density Standard 1.2471g/ml @25°C	100ml
DEN25190PY	Density Standard 1.3287g/ml @25°C	100ml
DEN25200PY	Density Standard 1.4117g/ml @25°C	100ml
DEN2515PY	Density Standard 1.5000g/ml @25°C	100ml
DEN25210PY	Density Standard 1.5775g/ml @25°C	100ml
DEN2516PY	Density Standard 1.6000g/ml @25°C	100ml
DEN25220PY	Density Standard 1.7441g/ml @25°C	100ml
DEN25230PY	Density Standard 1.9108g/ml @25°C	100ml
DEN25240PY	Density Standard 2.0775g/ml @25°C	100ml
DEN25250PY	Density Standard 2.2490g/ml @25°C	100ml
DEN25260PY	Density Standard 2.4175g/ml @25°C	100ml
DEN25270PY	Density Standard 2.5964g/ml @25°C	100ml
DEN25280PY	Density Standard 2.7493g/ml @25°C	100ml
DEN25290PY	Density Standard 2.9319g/ml @25°C	100ml
DEN25300PY	Density Standard 3.1043g/ml @25°C	100ml

Density Standards @ 40°C

Product No.	Description	Pack Size
DEN40010PY	Density Standard 0.6752g/ml @40°C	100ml
DEN40020PY	Density Standard 0.6872g/ml @40°C	100ml
DEN40030PY	Density Standard 0.6997g/ml @40°C	100ml
DEN40040PY	Density Standard 0.7109g/ml @40°C	100ml
DEN40050PY	Density Standard 0.7226g/ml @40°C	100ml
DEN40060PY	Density Standard 0.7343g/ml @40°C	100ml
DEN40070PY	Density Standard 0.7531g/ml @40°C	100ml
DEN40080PY	Density Standard 0.7733g/ml @40°C	100ml
DEN40090PY	Density Standard 0.7958g/ml @40°C	100ml
DEN40100PY	Density Standard 0.8207g/ml @40°C	100ml
DEN40110PY	Density Standard 0.8482g/ml @40°C	100ml
DEN40120PY	Density Standard 0.8945g/ml @40°C	100ml
DEN40130PY	Density Standard 0.9323g/ml @40°C	100ml
DEN40140PY	Density Standard 0.9857g/ml @40°C	100ml
DEN40150PY	Density Standard 1.0152g/ml @40°C	100ml
DEN40160PY	Density Standard 1.0642g/ml @40°C	100ml
DEN40170PY	Density Standard 1.1581g/ml @40°C	100ml
DEN40180PY	Density Standard 1.2408g/ml @40°C	100ml
DEN40190PY	Density Standard 1.3217g/ml @40°C	100ml
DEN40200PY	Density Standard 1.4039g/ml @40°C	100ml
DEN40210PY	Density Standard 1.5685g/ml @40°C	100ml
DEN40220PY	Density Standard 1.7339g/ml @40°C	100ml
DEN40230PY	Density Standard 1.8994g/ml @40°C	100ml
DEN40240PY	Density Standard 2.0649g/ml @40°C	100ml
DEN40250PY	Density Standard 2.2352g/ml @40°C	100ml
DEN40260PY	Density Standard 2.4028g/ml @40°C	100ml
DEN40270PY	Density Standard 2.5807g/ml @40°C	100ml
DEN40280PY	Density Standard 2.7329g/ml @40°C	100ml
DEN40290PY	Density Standard 2.9132g/ml @40°C	100ml
DEN40300PY	Density Standard 3.0852g/ml @40°C	100ml

Density Standards @ 50°C

Product No.	Description	Pack Size
DEN50010PY	Density Standard 0.6668g/ml @50°C	100ml
DEN50020PY	Density Standard 0.6791g/ml @50°C	100ml
DEN50030PY	Density Standard 0.6917g/ml @50°C	100ml
DEN50040PY	Density Standard 0.7033g/ml @50°C	100ml
DEN50050PY	Density Standard 0.7151g/ml @50°C	100ml
DEN50060PY	Density Standard 0.7269g/ml @50°C	100ml
DEN50070PY	Density Standard 0.7454g/ml @50°C	100ml
DEN50080PY	Density Standard 0.7653g/ml @50°C	100ml
DEN50090PY	Density Standard 0.7873g/ml @50°C	100ml
DEN50100PY	Density Standard 0.8118g/ml @50°C	100ml
DEN50110PY	Density Standard 0.8387g/ml @50°C	100ml
DEN50120PY	Density Standard 0.8868g/ml @50°C	100ml
DEN50130PY	Density Standard 0.9245g/ml @50°C	100ml
DEN50140PY	Density Standard 0.9777g/ml @50°C	100ml
DEN50150PY	Density Standard 1.0073g/ml @50°C	100ml
DEN50160PY	Density Standard 1.0562g/ml @50°C	100ml
DEN50170PY	Density Standard 1.1512g/ml @50°C	100ml
DEN50180PY	Density Standard 1.2346g/ml @50°C	100ml
DEN50190PY	Density Standard 1.3138g/ml @50°C	100ml
DEN50200PY	Density Standard 1.3973g/ml @50°C	100ml
DEN50210PY	Density Standard 1.5609g/ml @50°C	100ml
DEN50220PY	Density Standard 1.7257g/ml @50°C	100ml
DEN50230PY	Density Standard 1.8904g/ml @50°C	100ml
DEN50240PY	Density Standard 2.0551g/ml @50°C	100ml
DEN50250PY	Density Standard 2.2247g/ml @50°C	100ml
DEN50260PY	Density Standard 2.3916g/ml @50°C	100ml
DEN50270PY	Density Standard 2.5689g/ml @50°C	100ml
DEN50280PY	Density Standard 2.7207g/ml @50°C	100ml
DEN50290PY	Density Standard 2.9005g/ml @50°C	100ml
DEN50300PY	Density Standard 3.0721g/ml @50°C	100ml

Density Standards @ 60°C

Product No.	Description	Pack Size
DEN60010PY	Density Standard 0.6582g/ml @60°C	100ml
DEN60020PY	Density Standard 0.6708g/ml @60°C	100ml
DEN60030PY	Density Standard 0.6835g/ml @60°C	100ml
DEN60040PY	Density Standard 0.6955g/ml @60°C	100ml
DEN60050PY	Density Standard 0.7076g/ml @60°C	100ml
DEN60060PY	Density Standard 0.7196g/ml @60°C	100ml
DEN60070PY	Density Standard 0.7376g/ml @60°C	100ml
DEN60080PY	Density Standard 0.7572g/ml @60°C	100ml
DEN60090PY	Density Standard 0.7788g/ml @60°C	100ml
DEN60100PY	Density Standard 0.8027g/ml @60°C	100ml
DEN60110PY	Density Standard 0.8292g/ml @60°C	100ml
DEN60120PY	Density Standard 0.8790g/ml @60°C	100ml
DEN60130PY	Density Standard 0.9166g/ml @60°C	100ml
DEN60140PY	Density Standard 0.9695g/ml @60°C	100ml
DEN60150PY	Density Standard 0.9990g/ml @60°C	100ml
DEN60160PY	Density Standard 1.0478g/ml @60°C	100ml

Density Standards @ 80°C

Product No.	Description	Pack Size
DEN80010PY	Density Standard 0.6407g/ml @80°C	100ml
DEN80020PY	Density Standard 0.6538g/ml @80°C	100ml
DEN80030PY	Density Standard 0.6661g/ml @80°C	100ml
DEN80040PY	Density Standard 0.6798g/ml @80°C	100ml
DEN80050PY	Density Standard 0.6923g/ml @80°C	100ml
DEN80060PY	Density Standard 0.7047g/ml @80°C	100ml
DEN80070PY	Density Standard 0.7220g/ml @80°C	100ml
DEN80080PY	Density Standard 0.7407g/ml @80°C	100ml
DEN80090PY	Density Standard 0.7614g/ml @80°C	100ml
DEN80100PY	Density Standard 0.7844g/ml @80°C	100ml
DEN80110PY	Density Standard 0.8098g/ml @80°C	100ml
DEN80120PY	Density Standard 0.8629g/ml @80°C	100ml
DEN80130PY	Density Standard 0.9006g/ml @80°C	100ml
DEN80140PY	Density Standard 0.9520g/ml @80°C	100ml
DEN80150PY	Density Standard 0.9815g/ml @80°C	100ml
DEN80160PY	Density Standard 1.0302g/ml @80°C	100ml

Density Standards - Quality Range



Summary of Features & Benefits:

Commercial Benefits

- Extensive range (0.6538 - 1.0337 g/ml)
- No toxic heavy metals used in any formulation
- Can be used with any brand or type of vibrational density measuring instrument
- Presented in a high quality tamper proof amber glass bottle
- Customised formulations available

Technical Benefits

- Test results accredited to ISO17025. Uncertainty of measurement (assay procedure) $\pm 0.16\%$
- Produced in accordance with ASTM D4052 Guidelines
- Consistency of product - Independent, Traceable, Certified
- Certificates of Analysis and Safety Data Sheets available online
- ISO/IEC 17025 (INAB Ref:264T) Density Range 0.65-1.034g/ml

Reagecon manufactures an extensive range of Density Standards in accordance with ASTM D4052 for testing of Density, Relative Density and API Gravity of Liquids by Digital Density Meter. These materials can be used as calibration standards for density measurement by vibrational techniques or hydrometer based techniques.

The product range includes:

- 0.6960 - 1.0337g/ml @ 15°C
- 0.6619 - 1.0301g/ml @ 20°C
- 0.6878 - 1.0265g/ml @ 25°C
- 0.6752 - 1.0152g/ml @ 40°C
- 0.6668 - 1.0073g/ml @ 50°C
- 0.6582 - 0.9990g/ml @ 60°C
- 0.6538 - 1.0302g/ml @ 80°C



These products are prepared gravimetrically on a weight/weight basis. Both solute and solvent are weighed on a balance calibrated by Reagecon engineers using OIML traceable weights. Reagecon holds ISO 17025 accreditation for calibration of laboratory balances (INAB Ref:265C). The resulting Balance Certificate of Calibration is issued in accordance with the requirements of ISO/IEC 17025. The density of each standard is verified using a high performance calibrated density meter. The control of the density meter (identification no. - RRD015) is completed using high purity ISO Guide 34 accredited density standards similar in density value to these products.

Density Standards @ 15°C

Product No.	Description	Pack Size
DEN15010	Density Standard 0.6960g/ml @15°C	100ml
DEN15020	Density Standard 0.7073g/ml @15°C	100ml
DEN15030	Density Standard 0.7184g/ml @15°C	100ml
DEN15040	Density Standard 0.7298g/ml @15°C	100ml
DEN15050	Density Standard 0.7411g/ml @15°C	100ml
DEN15060	Density Standard 0.7524g/ml @15°C	100ml
DEN15070	Density Standard 0.7721g/ml @15°C	100ml
DEN15080	Density Standard 0.7933g/ml @15°C	100ml
DEN15090	Density Standard 0.8168g/ml @15°C	100ml
DEN15100	Density Standard 0.8428g/ml @15°C	100ml
DEN15110	Density Standard 0.8715g/ml @15°C	100ml
DEN15120	Density Standard 0.9135g/ml @15°C	100ml
DEN15130	Density Standard 0.9514g/ml @15°C	100ml
DEN15140	Density Standard 1.0040g/ml @15°C	100ml
DEN15150	Density Standard 1.0337g/ml @15°C	100ml

Density Standards @ 20°C

Product No.	Description	Pack Size
DEN20010	Density Standard 0.6919g/ml @20°C	100ml
DEN20020	Density Standard 0.7033g/ml @20°C	100ml
DEN20030	Density Standard 0.7148g/ml @20°C	100ml
DEN20040	Density Standard 0.7261g/ml @20°C	100ml
DEN20050	Density Standard 0.7374g/ml @20°C	100ml
DEN20060	Density Standard 0.7488g/ml @20°C	100ml
DEN20070	Density Standard 0.7683g/ml @20°C	100ml
DEN20080	Density Standard 0.7893g/ml @20°C	100ml
DEN20090	Density Standard 0.8126g/ml @20°C	100ml
DEN20100	Density Standard 0.8384g/ml @20°C	100ml
DEN20110	Density Standard 0.8668g/ml @20°C	100ml
DEN20120	Density Standard 0.9098g/ml @20°C	100ml
DEN20130	Density Standard 0.9476g/ml @20°C	100ml
DEN20140	Density Standard 1.0005g/ml @20°C	100ml
DEN20150	Density Standard 1.0301g/ml @20°C	100ml

Density Standards @ 25°C

Product No.	Description	Pack Size
DEN25010	Density Standard 0.6878g/ml @25°C	100ml
DEN25020	Density Standard 0.6993g/ml @25°C	100ml
DEN25030	Density Standard 0.7111g/ml @25°C	100ml
DEN25040	Density Standard 0.7223g/ml @25°C	100ml
DEN25050	Density Standard 0.7337g/ml @25°C	100ml
DEN25060	Density Standard 0.7452g/ml @25°C	100ml
DEN25070	Density Standard 0.7645g/ml @25°C	100ml
DEN25080	Density Standard 0.7853g/ml @25°C	100ml
DEN25090	Density Standard 0.8084g/ml @25°C	100ml
DEN25100	Density Standard 0.8340g/ml @25°C	100ml
DEN25110	Density Standard 0.8622g/ml @25°C	100ml
DEN25120	Density Standard 0.9060g/ml @25°C	100ml
DEN25130	Density Standard 0.9438g/ml @25°C	100ml
DEN25140	Density Standard 0.9969g/ml @25°C	100ml
DEN25150	Density Standard 1.0265g/ml @25°C	100ml

Density Standards @ 40°C

Product No.	Description	Pack Size
DEN40010	Density Standard 0.6752g/ml @40°C	100ml
DEN40020	Density Standard 0.6872g/ml @40°C	100ml
DEN40030	Density Standard 0.6997g/ml @40°C	100ml
DEN40040	Density Standard 0.7109g/ml @40°C	100ml
DEN40050	Density Standard 0.7226g/ml @40°C	100ml
DEN40060	Density Standard 0.7343g/ml @40°C	100ml
DEN40070	Density Standard 0.7531g/ml @40°C	100ml
DEN40080	Density Standard 0.7733g/ml @40°C	100ml
DEN40090	Density Standard 0.7958g/ml @40°C	100ml
DEN40100	Density Standard 0.8207g/ml @40°C	100ml
DEN40110	Density Standard 0.8482g/ml @40°C	100ml
DEN40120	Density Standard 0.8945g/ml @40°C	100ml
DEN40130	Density Standard 0.9323g/ml @40°C	100ml
DEN40140	Density Standard 0.9857g/ml @40°C	100ml
DEN40150	Density Standard 1.0152g/ml @40°C	100ml

Density Standards @ 50°C

Product No.	Description	Pack Size
DEN50010	Density Standard 0.6668g/ml @50°C	100ml
DEN50020	Density Standard 0.6791g/ml @50°C	100ml
DEN50030	Density Standard 0.6917g/ml @50°C	100ml
DEN50040	Density Standard 0.7033g/ml @50°C	100ml
DEN50050	Density Standard 0.7151g/ml @50°C	100ml
DEN50060	Density Standard 0.7269g/ml @50°C	100ml
DEN50070	Density Standard 0.7454g/ml @50°C	100ml
DEN50080	Density Standard 0.7653g/ml @50°C	100ml
DEN50090	Density Standard 0.7873g/ml @50°C	100ml
DEN50100	Density Standard 0.8118g/ml @50°C	100ml
DEN50110	Density Standard 0.8387g/ml @50°C	100ml
DEN50120	Density Standard 0.8868g/ml @50°C	100ml
DEN50130	Density Standard 0.9245g/ml @50°C	100ml
DEN50140	Density Standard 0.9777g/ml @50°C	100ml
DEN50150	Density Standard 1.0073g/ml @50°C	100ml

Density Standards @ 60°C

Product No.	Description	Pack Size
DEN60010	Density Standard 0.6582g/ml @60°C	100ml
DEN60020	Density Standard 0.6708g/ml @60°C	100ml
DEN60030	Density Standard 0.6835g/ml @60°C	100ml
DEN60040	Density Standard 0.6955g/ml @60°C	100ml
DEN60050	Density Standard 0.7076g/ml @60°C	100ml
DEN60060	Density Standard 0.7196g/ml @60°C	100ml
DEN60070	Density Standard 0.7376g/ml @60°C	100ml
DEN60080	Density Standard 0.7572g/ml @60°C	100ml
DEN60090	Density Standard 0.7788g/ml @60°C	100ml
DEN60100	Density Standard 0.8027g/ml @60°C	100ml
DEN60110	Density Standard 0.8292g/ml @60°C	100ml
DEN60120	Density Standard 0.8790g/ml @60°C	100ml
DEN60130	Density Standard 0.9166g/ml @60°C	100ml
DEN60140	Density Standard 0.9695g/ml @60°C	100ml
DEN60150	Density Standard 0.9990g/ml @60°C	100ml

Density Standards @ 80°C

Product No.	Description	Pack Size
DEN80020	Density Standard 0.6538g/ml @80°C	100ml
DEN80030	Density Standard 0.6661g/ml @80°C	100ml
DEN80040	Density Standard 0.6798g/ml @80°C	100ml
DEN80050	Density Standard 0.6923g/ml @80°C	100ml
DEN80060	Density Standard 0.7047g/ml @80°C	100ml
DEN80070	Density Standard 0.7220g/ml @80°C	100ml
DEN80080	Density Standard 0.7407g/ml @80°C	100ml
DEN80090	Density Standard 0.7614g/ml @80°C	100ml
DEN80100	Density Standard 0.7844g/ml @80°C	100ml
DEN80110	Density Standard 0.8098g/ml @80°C	100ml
DEN80120	Density Standard 0.8629g/ml @80°C	100ml
DEN80130	Density Standard 0.9006g/ml @80°C	100ml
DEN80140	Density Standard 0.9520g/ml @80°C	100ml
DEN80150	Density Standard 0.9815g/ml @80°C	100ml
DEN80160	Density Standard 1.0302g/ml @80°C	100ml

